

OPERATING AND SAFETY INSTRUCTIONS FOR



Set-Up

Plug the power cord from the adapter into the Mini-Mixer. Next plug the adapter into an outlet. The power switch for the control is located on the front of the unit. Once power is applied the green LED will light.

Loading the Mixer

Position up to 16 microcentrifuge tubes into the Mini-Mixer. The unit is now ready for operation.

Operating the Unit

Once the power switch is engaged, the mixer will start. Use the knob to adjust the frequency from 2 to 90 mixes per minute.

Maintenance

This mixer does not require regularly scheduled maintenance. However, regular inspection is recommended. Mixers should be protected from chemical spillage, mechanical damage and corrosive atmospheres so far as possible. Contamination, overheating and general misuse will greatly reduce the life of the Mini-Mixer.

Limited Warranty

We warrant products of manufacture to be free from defects in material and workmanship for one year and agree to repair or replace without charge any products

found defective upon examination at the factory. With proper care and operation, these products will give long and efficient service. Chemical spillage, overloading and general misuse will greatly reduce the service life. We are not responsible for damage to an apparatus due to improper installation or through attempts to operate the apparatus beyond its rated capacity, intentional or otherwise.

Limitations of Warranties

Apart from such written statement of warranty, there are no warranties, expressed, implied or statutory, which extend beyond the description of the products on the face hereof.

Specifications

- **Size:** 7-1/4"w x 3"d x 5-1/4"h
- **Electrical Rating:** (PIR) 100-240V, 50/60 Hz, 0.8A, (SEC) 6VDC, 2.0A with 6-ft, 2-wire cord
- **Weight:** 2.2 lbs.
- **Power Switch:** On/Off
- LED power indicator
- **Frequency of mix:** Varies from 2 to 90 mixes per minute
- **Ambient Temperature:** 77°F/25°C +/-5°
- **Operating Environment:** 0 to 90% relative humidity 32 to 100°F/0 to 37°C



MTC • MicroSolv Technology Corporation

1 Industrial Way West, Bldg E, Eatontown, NJ 07724 USA

Ph. 1-732-578-1777 • Fx. 1-732-578-9777

www.mtc-usa.com Email: info@mtc-usa.com