



## To obtain greater sensitivity than Na adducts for sugars could you add ammonia NH<sub>4</sub> instead - FAQ

In LCMS, you will get much better sensitivity of Na-sugar adducts in the EICs compared to sugars themselves.

The same principle will work for NH<sub>4</sub> adducts as well, but there are notable drawbacks. The problems with NH<sub>4</sub> is that it will change the pH of the mobile phase and its concentration is hard to control in the sample. Hence, given both strategies will work, there is no reason to choose ammonia NH<sub>4</sub> over sodium Na.



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)