



Equilibrating HPLC Column Suggestions to Protect the Columns - Tips & Suggestions

Date: 22-APRIL-2012 Last Updated: 1-DECEMBER-2025

Always ensure that the column is fully equilibrated with your starting mobile phase before making any injections.

When you are equilibrating your columns for the first time or whenever changing your HPLC methods it is wise to consider not putting anything into the columns that can precipitate or is immiscible with the storage or current solvent in the column.

Depending on column type, equilibrating could range from 10 to 15 column bed volumes for typical reversed phase to 50 to 100 bed volumes with polar embedded phases or HILIC columns.

Beware of buffers that might be in your column before you change the mobile phase composition with respect to the organic content. This can cause immediate precipitation and ruin your HPLC columns.



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com