

## Low pH buffers Used to Separate Acidic Compounds in RP - HPLC Primer

*Date: 22-APRIL-2012 Last Updated: 1-DECEMBER-2025*

By suppressing their ionization at low pH, acids will be more like neutral compounds and will retain on Reversed Phase HPLC columns much the same way that neutral compounds will. The same is true for using bases with basic compounds.

Another benefit of running on Type A and Type B silica at low pH is that the silanol activity is reduced. This should sharpen your peaks.



Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)