

Using the wavelength calibration solution holmium oxide for HPLC UV detector performance qualification - Tips & Suggestions

Date: 14-APRIL-2012 Last Updated: 2-JULY-2025

Performing the wavelength accuracy procedure from our HSQ™ and PQ™ Qualification Kits correctly is very easy and requires very little qualification solution.

While pulling a vacuum and filling the flow cell it is possible that you may see small bubbles coming through the plastic tubing. These bubbles seem to come from the tubing connections. Pulling solution through over and over again, waiting for them to disappear will use up time and the holmium oxide (HOX) solution. Be sure to have good solid tubing connections and pull the solution through the cell *slowly and evenly*. The bubbles will not dissipate on their own accord.

One just has to make sure there are no bubbles in the cell; so just pull a drop through, break the vacuum then take a spectrum. If by chance there is a small bubble in the cell, and you get a bad spectrum then simply pull another drop through and try again.

For example: One drop of holmium oxide basically should be enough to flush a 10µL cell completely with the solution. It is very rare that an experienced user does not obtain a good quality spectrum the first time.

Click [HERE](#) for Chemical Solutions DIY Qualification Kits™ Ordering Information and Pictures.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com