



Since Cogent TYPE-C columns can be used in Reversed Phase or Normal Phase what do you suggest to switch these solvent systems which are not miscible - FAQ

**Suggested Procedure:**

**A** – Changing from Reversed Phase to Normal Phase HPLC; pump 100% methanol for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methylene chloride. The column is ready to be equilibrated with mobile phase for NP HPLC.

**B** – Changing from Normal Phase to Reverses Phase HPLC; pump 100% methylene chloride for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methanol. The column is ready to be equilibrated with mobile phase for RP-HPLC.

A simple 30 minute procedure allows switching from one mode to another.



Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)