



Ammonium Acetate in LCMS Requires Attention to Cleaning - Tips & Suggestions

Date: 14-APRIL-2020 Last Updated: 1-DECEMBER-2025

Ammonium Acetate is a wonderful additive for LCMS using HILIC or Aqueous Normal Phase (ANP) methods. However, be aware that even in low concentrations, the inlet source of the MS will become clogged with over time and will require cleaning.

If you notice that your peaks are disappearing or the integration of your desired peaks decrease, it is time to perform some preventative maintenance.

[Cogent Diamond Hydride Product Page](#)



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com