

Dimensions: 2.1 x 150 mm

Mobile Phase:

50% A: DI Water / 0.1% Formic Acid

50% B: Acetonitrile / 0.1% Formic Acid

Injection vol.: 1 µL

Flow rate: 0.5 mL / minute

Detection: ESI – POS - PerkinElmer, Flexar SQ 300 Mass Spectrometer

Temperature: 25°C

Sample Preparation:

- Standard: 5 ppm of Limonin in 20% DI Water / 0.1% Formic Acid / 40% Acetonitrile / 40% Methanol.
- Spiked Orange Juice preparation: Orange juice was spiked with 2.5 ppm Limonin, filtered, and injected.

t_o : 1.1 minute

Note: Limonin is a bitter compound which may negatively affect juice quality. The compound is found in the seeds and membrane tissue of the fruit. It is very important for groves to determine the level of Limonin in juice so the correct recovery settings for the juice production can be set. The level of Limonin can change dramatically from season to season. It also depends on the fruit size. The analysis of Limonin is crucial in production of high quality non bitter fruit juices.



Attachment No 281 Limonin in Orange Juice LCMS pdf 0.2 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com