

Standard Silicone pre Slit Sealing Mat Compared to Solvent Resistant Mat - Tips & Suggestions

Date: 7-NOVEMBER-2020 Last Updated: 27-OCTOBER-2025

Comparison of 96 well sealing mats, Silicone Rubber vs. Chemical Resistant U-2D mats.

Silicone rubber mats should only be used with pure aqueous and buffer solutions.

The solvent resistant mat can be used with methanol, acetonitrile and water / buffers. These should not be used with non-polar solvents or aggressive acid or base such as ethyl acetate. The maximum temperature for these mats is 100°C.

[U-2D Silicone Mats Product Page](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com