

What is the difference between a Cap Liner and a Cap Septum - HPLC Primer

21-MAY-2021 Last Updated: 20-JUNE-2025


In chromatography and laboratory vial systems, cap liners and cap septa serve different purposes, even though they may appear similar in material or thickness.

Cap Liner

- A liner is typically installed in a solid-top cap (i.e., a cap without a hole).
- It cannot be pierced and is not designed for needle access.
- Its primary function is to seal the vial, preventing sample evaporation or contamination during storage or transport.
- Liners may be made from materials such as PTFE, silicone, or polyethylene, and are selected based on chemical compatibility.

Cap Septum

- A septum (plural: septa) is installed in an open-top cap, which has a hole to allow needle access.
- It is designed to be pierced by an autosampler needle or syringe during sample injection.
- After piercing, the septum reseals itself, maintaining vial integrity and minimizing exposure to air or contaminants.
- Septa are often made from layered materials (e.g., PTFE/silicone) to balance chemical resistance and resealability.

 **Fun Fact:** The term *septum* comes from Latin, meaning “a wall” or “partition.” In chromatography, it acts as a barrier that protects the sample until it's accessed.

Why It Matters

- Using the correct closure type is essential for sample integrity, instrument compatibility, and analytical accuracy.
- Always match the cap type (solid or open-top) with the appropriate liner or septum based on your application—whether for storage or injection.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com