


How to Complete the BASIK™ Mobile Phase Filter Degasser Assembly - How To

Date: 18-JULY-2025 Last Updated: 30-NOVEMBER-2025

To complete the BASIK™ Mobile Phase Filter Degasser assembly:

- The **glass membrane support** (Catalog No. [58800-47-GS](#)) is the base component that holds the filter membrane in place.
- A **47 mm membrane filter disc** is placed **on top of this support**. (Via Red Circle, below.)

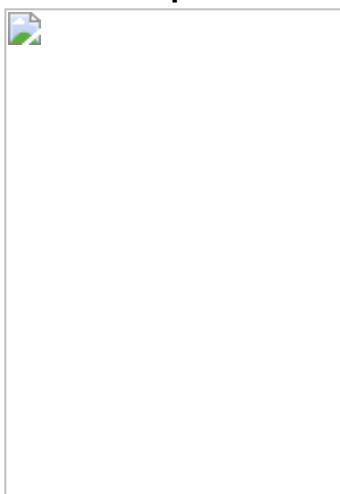

- The **glass funnel** (Catalog No. [58800-47-GF](#)) is then positioned **above the filter**, allowing solvent to pass through the membrane during filtration.
- To secure the funnel and membrane support together, use the **blue metal clamp** (Catalog No. [58800-47-CL](#)).



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com