


Ciprofloxacin Analyzed With LCMS - AppNote


Retention with Good Peak Shape for this Polar Compound

Ciprofloxacin has several amine groups and therefore is susceptible to peak tailing in Reversed Phase methods. The USP assay method for Ciprofloxacin tablets uses Triethylamine as an Ion Pair Agent to reduce tailing and in addition, Phosphoric Acid is included in the Mobile Phase.

These conditions make the USP method incompatible with MS detection and therefore limit its applications. In this method, only Formic Acid is required in order to obtain a peak that meets the system suitability specification for the API tailing factor. This Method demonstrates how various Type-C™ application notes can be successfully adapted for LCMS.

The Peak Shape for this polar compound meets the System Suitability specification for the USP Tailing Factor.

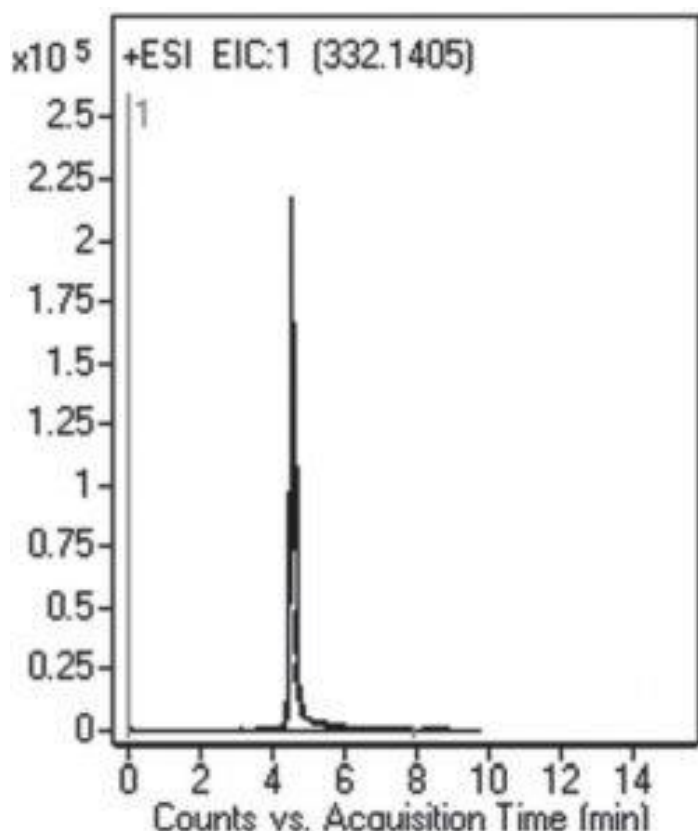

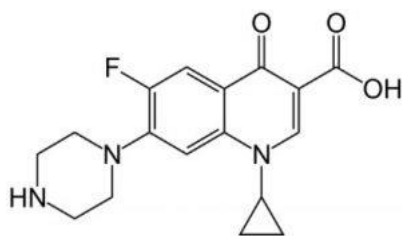

Peak:

Ciprofloxacin 332.1405 m/z (M+H)⁺

Method Conditions

Column: Cogent Diamond Hydride™, 4μm, 100Å

Catalog No.: [70000-15P-2](#)

Dimensions: 2.1 x 150mm

Mobile Phase:

A: 50% DI Water / 50% 2-Propanol / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

Gradient:

Time (minutes)	%B
0	90
3	10
6	10
7	90

Post Time: 3 minutes

Injection vol.: 1μL

Flow rate: 0.4mL / minute

Detection: ESI – pos – Agilent 6210 MSD TOF Mass Spectrometer

Sample Preparation: A 250mg strength Ciprofloxacin tablet was ground and dissolved in 25mL of 50:50 Solvent A / Solvent B diluent. Solution was sonicated 10 minutes and filtered through 0.45μm Nylon Syringe Filter (MICROSOLV Tech Corp.). Sample for injection was diluted 1:100 with the same diluent.

t₀: 0.9 minutes

Note: Ciprofloxacin is a Fluoroquinolone antibiotic which acts by inhibition of DNA Gyrase and Topoisomerase IV. It was once used as a drug of last resort for bacterial infections, but accumulated resistance has made the drug less effective in recent years.



Attachment

No 182 Ciprofloxacin Analyzed with LCMS pdf 0.2 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com