

Evaporation Study Soft-Guard Caps - Tech Information

Date: 15-JANUARY-2019 Last Updated: 2-FEBRUARY-2026

Evaporation Study: Soft-Guard™ vs. Market-Leading Pre-Slit Caps

Purpose: Compare evaporation rates between Soft-Guard™ caps and market-leading / private-label caps with pre-slit septa.

1. Materials & Methods

- **Sample Setup:**
 - 10 vials labeled #1–#10
 - 5 Market Leader (M) caps and 5 Soft-Guard™ (SG) caps
 - AQ Brand Maximum Recovery screw-top vials used
 - **Procedure:**
 - Empty vials with caps weighed (tare weight, w_{tare})
 - 800 μ L methanol added; vials weighed again (w_0)
 - Vials stored at room temperature for ~161 hours
 - Final weight recorded (w_{161})
-

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
2. Visual Observations

- **Market-Leading Caps:**
 - Septa appeared flat when loose
 - Became concave when attached to vials
 - **Soft-Guard™ Caps:**
 - Loose caps showed convex septa with slightly open slit (“smiling caps”)
 - When tightened, slit closed and surface appeared flat to slightly convex
-

3. Results

- Market-leading pre-slit caps lost **twice as much solvent** compared to Soft-Guard™ caps, despite Soft-Guard™ caps initially showing more exposed slits when loose.

Cap Type	Average % Loss
Market Leader (M)	5.4%
Soft-Guard™ (SG)	2.4%



Key Takeaways : Soft-Guard™ caps significantly reduce evaporation compared to market-leading alternatives. Design features (slit closure upon tightening) contribute to improved performance.



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com