
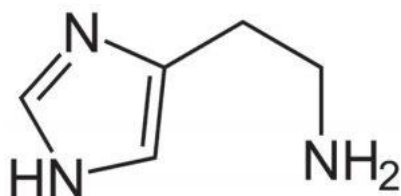


Histamine in red and white wine analyzed with LCMS - AppNote

LCMS Method without Derivatization

This Method is simple, fast, robust, and does not require derivatization, which is the main advantage over other HPLC Methods used in the Assay of this polar compound. It was demonstrated that Histamine is not present in the White Wine sample (*Figure B*). Sample preparation was minimal and involved just filtration using syringe filters. The sample of Red Wine contained Histamine at a level that was determined to be below 10mg / L (*Figure C*).

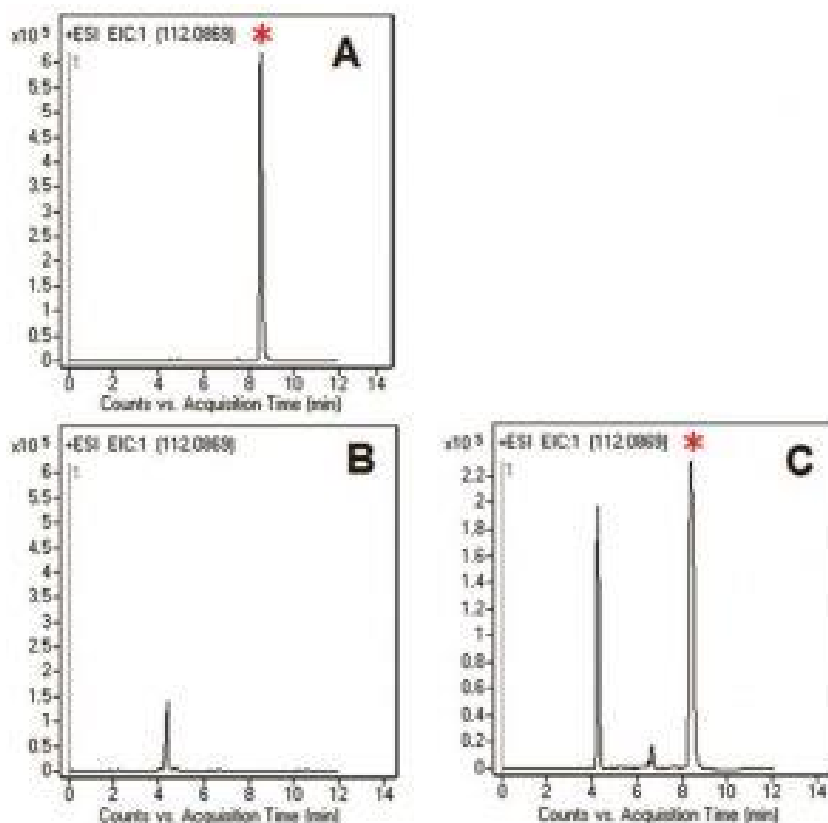

Method Conditions

Column: Cogent Diamond Hydride™, 4μm, 100Å

Catalog No.: [70000-15P-2](#)

Dimensions: 2.1 x 150mm

Mobile Phase:

A: 50% DI Water / 50% 2-Propanol / 0.1% Formic Acid

B: Acetonitrile / 0.1% Formic Acid

Gradient:

Time (minutes)	%B
0	80
5	10
7	10
8	80

Post Time: 5 minutes (can be reduced for standards)

Injection vol.: 1 µL

Flow rate: 0.4 mL / minute

Detection: ESI – POS - Agilent 6210 MSD TOF Mass Spectrometer

Figures:

A: *Histamine Standard*. 1 mg / mL Histamine Dihydrochloride in DI Water / Methanol 80:20. The solution was filtered through a 0.45 µm Filter (MICROSOLV Tech Corp.). Sample for injection was diluted 1:100 with 50:50 Solvent A:B mixture. * = **Histamine Peak**

B: *White wine*. The wine was filtered through a disposable 0.45 µm Filter (MICROSOLV Tech Corp.) and injected as is. No histamine peak. Extra peak is from wine matrix.

C: *Old red wine*. The wine was filtered through a disposable 0.45 µm Filter (MICROSOLV Tech Corp.). Sample for injection was diluted 1:5 with 50:50 Solvent A:B mixture. * = **Histamine Peak**
Extra peaks are from wine matrix.

t₀: 0.9 minutes

Note: Histamine is known for its role in allergic response. Release of Histamine plays a role in inflammation, gastric acid secretion, microcirculation and neurotransmission in mammalian brains. Measurement of Histamine levels in body fluids has been used in clinical analysis in various diseases such as Preeclampsia, Asthma, Cancer, Mastocytosis and in the progression of Periodontitis. Histamine is also present in many foods and beverages, especially red wine and spoiled food. Ingesting histamine can cause migraines, sweating and nausea.



Attachment

Histamine in red and white wine analyzed with LCMS pdf [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com