



Coricidin HBP Day Tablet Analyzed by HPLC- AppNote

Formulation of Guaifenesin and Dextromethorphan HBr

Click [HERE](#) for Column Ordering Information.

This Method shows the separation of two active ingredients in a formulation of Guaifenesin and Dextromethorphan. Dextromethorphan often exhibits tailing in many HPLC methods, but here an excellent Peak shape is observed.

This Method is suitable for routine assay of these compounds in combination formulations. A change in wavelength was used in order to obtain similar Peak heights between the two compounds, since their concentrations differ significantly.

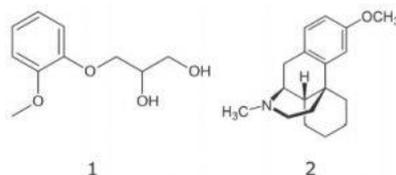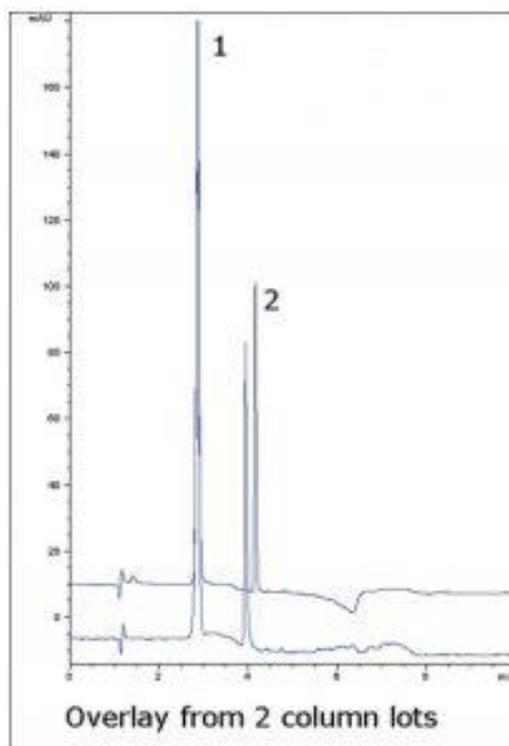

Peaks:

1. Guaifenesin, 2. Dextromethorphan

Method Conditions

Column: Cogent Phenyl Hydride™, 4 µm, 100Å

Catalog No.: 69020-75P

Dimensions: 4.6 x 75 mm

Mobile Phase:

A: DI Water / 0.1% TFA (v/v)

B: Acetonitrile / 0.1% TFA (v/v)

Gradient:

Time (minutes)	%B
0	20
1	20
5	90
6	20

Temperature: 35°C**Post Time:** 4 minutes**Flow rate:** 1.0 mL / minute**Detection:** UV @ 254 nm (0-3.5 minutes), 215 nm (3.5-10 minutes)**Injection vol.:** 2 µL

Sample Preparation: Coricidin High Blood Pressure (HBP) Day® tablet containing 200 mg Guaifenesin and 10mg Dextromethorphan HBr was ground and added to a 50mL volumetric flask. A portion of 50:50 Solvent A / Solvent B was added and the flask was sonicated 10 minutes. Then it was diluted to mark and mixed well. A portion was filtered with a 0.45 µm Nylon Syringe Filter (MICROSOLV Tech Corp.). This filtrate was used for injections. The Peak identities were confirmed with individual standards.

t_o: 0.9 minutes

Note: *Guaifenesin is an expectorant agent and Dextromethorphan is an antitussive. This particular cold and cough formulation is designed for individuals with high blood pressure.*



Attachment No 206 Coricidin HBP Day Tablet Analyzed by HPLC pdf 0.3 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.comWebsite: www.mtc-usa.com