
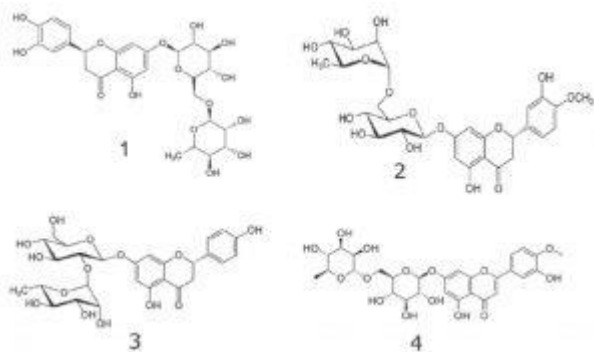
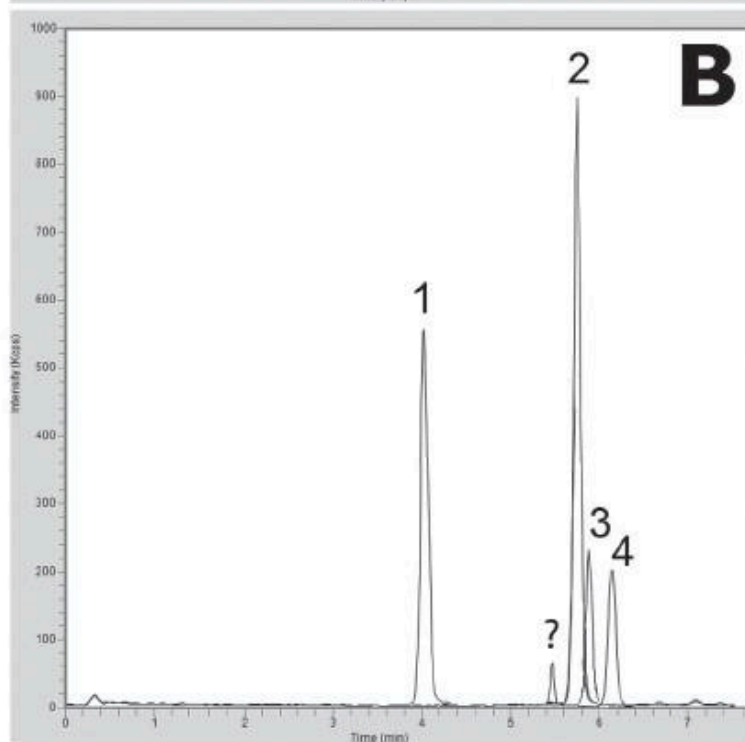


Spiked Hesperidin Extract Analyzed with LCMS - AppNote

Eriocitrin, Hesperidin, Naringin, Diosmin

Click [HERE](#) for Column Ordering Information.

In this method, the Hesperidin extract was spiked with three glycosidic flavonoids. The Cogent Phenyl Hydride Column was used to separate the main component of the extract (Hesperidin) and three standards. The MS detection helped to distinguish between the Naringin peak of the standard and a small impurity peak (peak marked "?"). When the Hesperidin extract sample was analyzed, only peaks for Hesperidin and a small peak for Diosmin were detected.

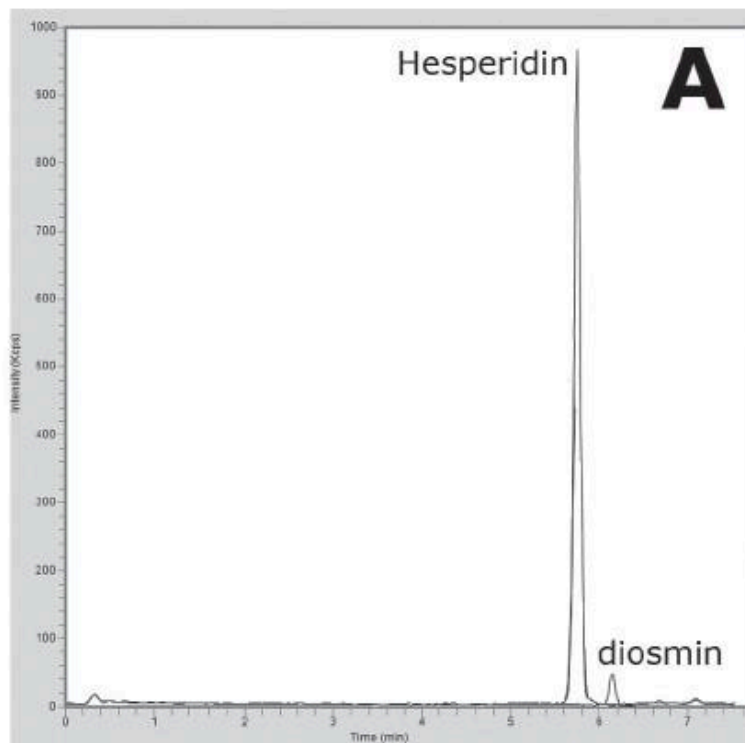

Peaks:

1. Eriocitrin 595.2 [M+H]⁺

2. Hesperidin 541.2 m/z [M+H]⁺
3. Naringin 581.2 m/z [M+H]⁺
4. Diosmin 609.2 m/z [M+H]⁺

Method Conditions

Column: Cogent Phenyl Hydride, 4µm, 1 00Å

Catalog No.: [69020-05P-2](#)

Dimensions: 2.1 x 50 mm

Mobile Phase:

A: DI Water / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

Gradient:

Time (minutes)	%B
0	10
1	10
9	70
10	10

Post Time: 3 minutes

Injection vol.: 1µL

Flow rate: 0.4 mL/minute

Detection: ESI – POS - PerkinElmer, Flexar SQ 300 Mass Spectrometer

Samples:

A: Hesperidin extract (1mg/mL) in DI Water, diluted 1:100 with Solvent B

B: Hesperidin extract (1mg/mL) in DI Water, diluted 1:100 with Solvent B and spiked with 20 ppm of the standards (Eriocitrin, Naringin, Diosmin)

Note: *Flavonoids are phenolic substances isolated from a variety of plants and fruits. They are low molecular weight compounds and in plants they are usually bound to sugar molecules as O-glycosides. Flavonoids are present in foods and beverages of plant origin, such as fruits, vegetables, tea, wine etc. They can be divided into fourteen classes, including flavonols, flavones, anthocyanins, isoflavonoids, and flavanones. They have been shown to have anti-inflammatory, anti-allergic, anti-carcinogenic, antihypertensive and anti-arthritis, and free radical scavenging properties.*



Attachment

No 293 Spiked Hesperidin Extract pdf 0.3 Mb [Download File](#)

Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com