


## Nucleoside Prodrug Analogs Analyzed with HPLC - AppNote

### Polar and Non-Polar Anticancer Prodrugs Analyzed in the Same Run

In this Method, we display Retention of both polar and nonpolar compounds in a simple Isocratic run. Adjustment of the aqueous to organic ratio can be used to make either the polar or more hydrophobic compounds elute first.

In the example presented, the more hydrophobic compounds are retained longer but retention and separation of the polar constituents is still achieved. If greater retention of the polar compounds is desired then a higher percentage of Acetonitrile is required.

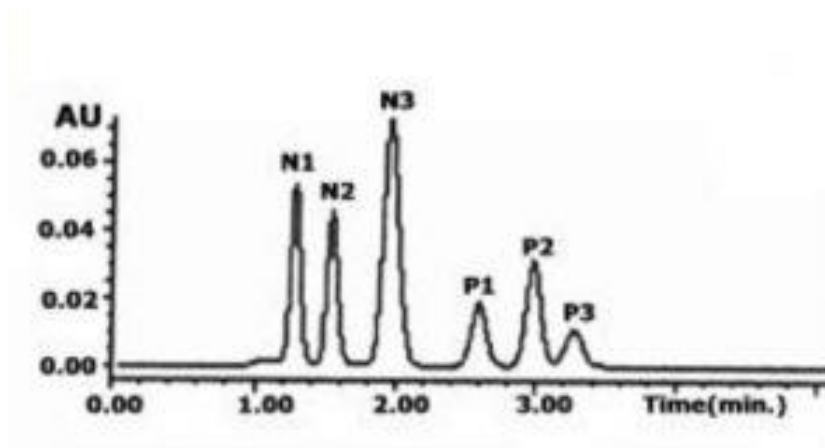

### Method Conditions

**Column:** Cogent UDC Cholesterol™, 4µm, 100Å

**Catalog No.:** 69069-7.5P

**Dimensions:** 4.6 x 75mm

**Mobile Phase:** 60% Acetonitrile / 40% DI Water / 0.5% Formic Acid

**Injection vol.:** 5µL

**Flow rate:** 1 mL / minute

**Detection:** UV @ 270nm

**Sample Preparation:** Proprietary Drugs. N1-N3 Nucleoside Analogues. P1-P3 have greater hydrophobicity with polar functional groups. 1µg / mL of each dissolved in the Mobile Phase

---

**Notes:** Nucleoside analogue anticancer “pro” drugs are difficult to retain on conventional HPLC Columns because of their highly polar natures. As such, Ion Pair Reagents are used in order to achieve sufficient Retention. These additives interfere with MS detection and are not preferred. With this Method, you can achieve good Retention paired with a MS-Compatible Mobile Phase.

---



**Attachment No 05 Nucleoside Analogs Analyzed with HPLC pdf** 0.1 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)