
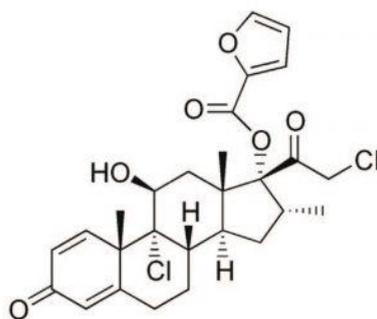


## Mometasone Furoate in Topical Cream Analyzed with HPLC - AppNote

### Separation of Prodrug from Acid Hydrolysis Products

This Method shows the Separation of an ester Prodrug, Mometasone Furoate, from another compound formed during acid hydrolysis. This Peak is quite possibly the active form, Mometasone. The two Chromatograms show the non-degraded extract (*Figure A*) and an acid hydrolysis extract (*Figure B*).

Hydrolysis of the ester group would make the resulting Mometasone API less hydrophobic, and therefore earlier elution is predicted. Indeed, this is what is observed in the data. An excipient Peak from the cream matrix is also observed just after the solvent front and does not interfere with either the Prodrug peak or the API.

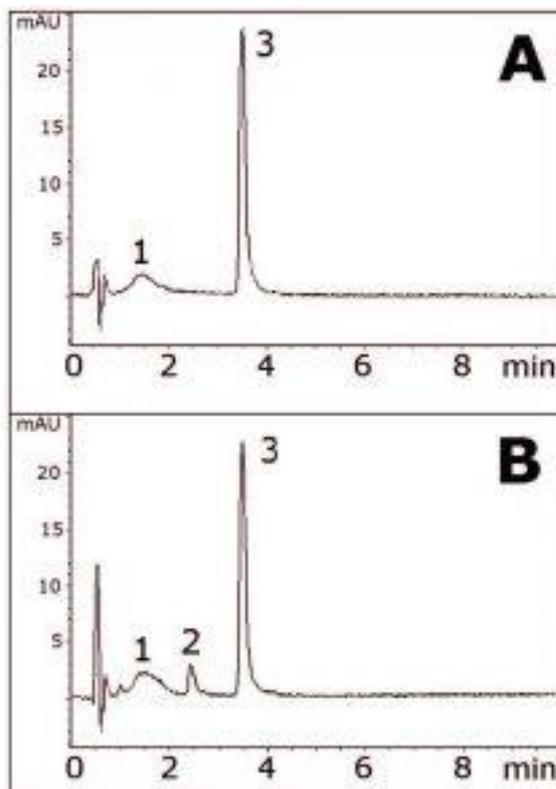

Mometasone Furoate

## Peaks:

1. Excipient,
2. Degradant (possibly Mometasone),
3. Mometasone Furoate (Prodrug)

## Method Conditions

**Column:** Cogent Bidentate C8 2.0™, 2.2 µm, 120 Å

**Catalog No.:** [40208-05P-2](#)

**Dimensions:** 2.1 x 50 mm

**Mobile Phase:** 50:50 DI Water / Acetonitrile (v/v)

**Injection vol.:** 2 µL

**Flow rate:** 0.3 mL / minute

**Detection:** UV @ 248 nm

**Sample Preparation:** 5.0 g of a 0.1% Mometasone Furoate USP topical cream was weighed in a 250 mL Erlenmeyer flask with a stirring bar. 50 mL Methanol was added and the flask was stirred for 1 hour with a stopper covering the flask. Then a portion was filtered with a 0.45 µm Nylon Syringe Filter (MICROSOLV Tech Corp.). Next, one portion of the filtrate was diluted 1:5 with Methanol (*Figure A*) and a second portion was diluted 1:5 with a diluent of 90:10 Methanol / 1N HCL and heated in a dry bath at 80°C for 10 minutes. (*Figure B*).

---

*Note:* Mometasone Furoate is a prodrug where the API is formed by hydrolysis of the ester group. It is used to reduce inflammation due a variety of ailments and is available under trade names such as Elocon®, Novasone®, and Asmanex®. In addition to a topical ointment, the drug is also available as a dry powder inhaler or a nasal spray depending on the application.

---



**Attachment No 325 Mometasone Furoate in Topical Cream Analysis pdf 0.3 Mb** [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)