

Identify Used U-2D Inserts for Multiple Projects or in Small Sample Sets - Tips & Suggestions

You can distinguish U-2D™ inserts (used vs. unused) by the following ways:

- A. Used samples can be indicated on the silicone mat/film using a marker.
- B. Alternatively, the film or mat can be cut with scissors to an appropriate size to protect unused inserts or to help keep track of used inserts.
- C. Users can selectively remove unused (or used) inserts as needed. Inserts should be carefully handled/stored to prevent contamination prior to use or transfer.

Click [HERE](#) for U-2D™ Micro-Sample Management System Ordering Information

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com