

## OBSOLETE - What maximum flow rate can I use for a PTFE 10m 0.75mm ID Knitted Reactor Coil KRC with isopropanol - FAQ

**Question:** I am using a PTFE Knitted Reactor Coil KRC with 10m length and 0.75mm ID with isopropanol as the fluid I am pumping. What is the maximum flow rate I can use under these conditions?

**Answer:** At ambient temperature, you should operate at no more than 10mL / minute with isopropanol in order to stay within the mechanical stability specifications of the KRC tubing. Ramp up the flow rate slowly from zero in order to avoid a sudden high pressure.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)