



Cogent Column Filters Do Not Need Special Fittings to use with HPLC column - Tech Information

Date: 22-APRIL-2012 Last Updated: 3-NOVEMBER-2025

The pre-column, Cogent Column Filter is completely compatible with all 1/16th", 10-32 internal fitting ports . Most modern brands of HPLC columns are fully compatible.

Finger tighten the filters right into the inlet side of the guard or analytical column (universal fit) and a finger tight male nut to hold your stainless or PEEK tubing inserted into the inlet side of the filter. Add 1/8 of a turn once you feel the filter is fairly tightly installed for the best seal.

[Click here to view more information about Column Protection](#)



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com