

What is the Maximum Liquid Volume for a 25mm AQ Brand Syringe Filter - FAQ

Date: 22-APRIL-2012 Last Updated: 3-NOVEMBER-2025

You can use up to 100ml and sometimes more, in a single 25mm AQ™ syringe filter if you are using an aqueous solvent and your sample is not cloudy.

However, the best way to determine if a filter should be replaced from service is when you get a very fast drop in back pressure or a constant rise in back pressure. Both indicate it is time to replace them. Reusing the same syringe filters for multiple samples is never recommended due to carry over problems.

Click [HERE](#) for syringe filters ordering information and pictures.

Attachments

MICROSOLV filters equivalency study pdf [Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com