

What Type of Fitting is Best to Use with the HPLC and LCMS Columns - FAQ

Date: 18-FEBRUARY-2015 Last Updated: 3-NOVEMBER-2025

We generally recommend the Direct Adaptive Di-Ad™ column connectors instead of a fitting that can be re-swaged each time one of our columns is removed (*both standard 4um and 2.0™ phases*) to enable a clean connection when changing HPLC columns. The Di-Ad™ allows proper and easy installation of Cogent columns as well as different columns without the hassles of swaged ferrules.

We do not recommend Di-Ad™ or PEEK fittings for UHPLC.

Click [HERE](#) for Di-Ad™ column connector information and pictures.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com