



Is the Cogent Guard Column Holder Supposed to Close All The Way - FAQ

Date: 6-JULY-2015 Last Updated: 30-NOVEMBER-2025

When you insert the guard column cartridge and close the holder, you may notice that the holder does not appear to close all the way.

This is normal and within the design specifications of the unit. With the cartridge placed in the holder, simply give the holder a quarter turn with a wrench and this should suffice to provide adequate sealing of the guard column.



Printed from the Chrom Resource Center
Copyright 2025, All Rights Apply
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900
Fax: (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com