



## Closing the Cogent Guard Column Holder - FAQ

*Date: 6-JULY-2015 Last Updated: 30-NOVEMBER-2025*

When you insert the guard column cartridge and close the holder, you may notice that the holder does not appear to close all the way.

This is normal and within the design specifications of the unit. With the cartridge placed in the holder, simply give the holder a quarter turn with a wrench and this should suffice to provide adequate sealing of the guard column.



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)