

Blockage Minimization in HPLC Columns - HPLC Primer

Date: 5-NOVEMBER-2015 Last Updated: 23-FEBRUARY-2026

Overview

Blockages in HPLC columns can lead to high backpressure, distorted peak shapes, and compromised analytical performance. Preventing these issues is far easier and more effective than attempting to resolve them after they occur.

Understanding the common sources of particulate and chemical contamination allows users to implement strategies that significantly extend column life and maintain method reliability.

1. Sample-Related Causes and Prevention

Complex samples—such as plasma, biological fluids, food extracts, or environmental materials—often contain particulate debris and matrix components that can accumulate on the column inlet frit.

Key Prevention Steps

- Use a 0.45 μm or 0.2 μm syringe filter before transferring sample into autosampler vials to prevent particulate contamination.
 - Employ Cogent Column Filters to stop particulates from reaching the column frit and causing irreversible blockage.
 - Use a guard column to retain matrix components that may chemically interact with the stationary phase, preserving column performance and longevity.
-

2. Mobile Phase Cleanliness

The mobile phase itself can introduce contamination, especially when buffers, salts, or microbial growth are present.

Best Practices

- Vacuum-filter all mobile phases before pouring them into solvent reservoir bottles to remove dust, bacteria, undissolved buffer crystals, and particulate matter.
 - Ensure buffer salts are fully dissolved before filtration and use proper solvent preparation procedures to minimize precipitation.
 - Replace mobile phase regularly and store aqueous phases in clean, bacteria-free conditions.
-

3. Instrument-Generated Contaminants

Particulates do not always originate in the sample—they can also be generated by the HPLC instrument itself.

Common Sources

- Wear from pump seals
- Injector rotor debris
- Valve particulates
- Aging internal tubing shedding material

Prevention Tip

Install a Cogent Column Filter to capture these internal particulates before they reach the column, providing ongoing protection against instrument-generated contamination.

Ordering Information

Click the links below for product details and images:

- Cogent Column Filters – Click [HERE](#) for ordering information & pictures
- Guard Columns – Click [HERE](#) for ordering information & pictures
- AQ Brand Syringe Filters – Click [HERE](#) for ordering information & pictures



Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com