

## How does the shorter column supplied in the HSQ kit affect the data compared to the longer column from the PQ Kit - FAQ

*Date: 9-APRIL-2014 Last Updated: 1-JULY-2025*

**The HSQ Kit™ comes with a 4.6 x 75mm column that is used in the qualification procedure.**

The older version of the the HSQ Kit™ (called the PQ Kit™) comes with a 4.6 x 150mm column instead. This difference will not affect the qualification data results and the columns can be used **interchangeably**. The only difference for the end-user is that the HPLC runs take much less time due to the shorter column length when using the HSQ Kit™.

Hence, this a great advantage over the older version.

Click [HERE](#) for HSQ & PQ kits ordering information and pictures.

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)