



Different Tests in Chemical Solutions HSQ Kit Uses the Same Method Conditions - Tech Information

Date: 14-MAY-2014 Last Updated: 28-JUNE-2025

All tests included in both of the Chemical Solutions HSQ™ Qualification Kits for HPLC are designed to run under the same method conditions. This consistency allows for seamless overnight operation using an autosampler.

To perform Performance Qualification (PQ) validation, simply configure your system with the following parameters and let it run:

Wavelength:	273 nm
Injection Volume:	10 µl (except for volume linearity)
Flow Rate:	1 mL / minute
Temperature:	Ambient

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com