



## Metformin Impurities Melamine and Cyanoguanidine Determination in HCL Formulation - AppNote


### Simple Separation of API from Melamine and Cyanoguanidine

A simple Method was developed for the analysis of the widely prescribed anti-diabetic drug Metformin Hydrochloride and for determination of two impurities, Cyanoguanidine and Melamine, in tablet formulations.

In addition, to Accurate Mass all three compounds, the Peaks were confirmed by injections of standards. The USP specified impurity limits are not more than 0.02% and 0.01% for Cyanoguanidine and Melamine respectively.

The precision of this Method was evaluated by calculating %RSD of the Peak areas of five injections (see *Figure B*). The obtained value was 0.2%.

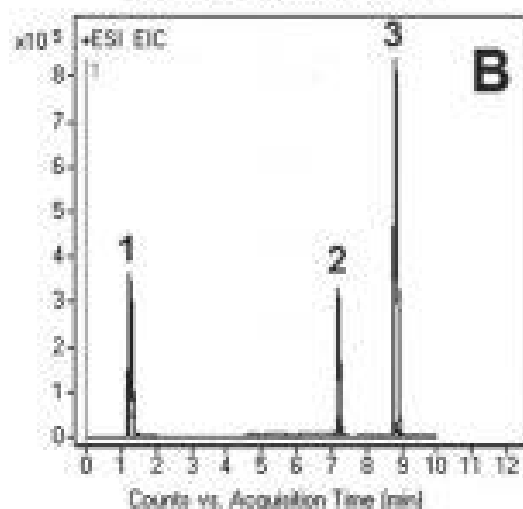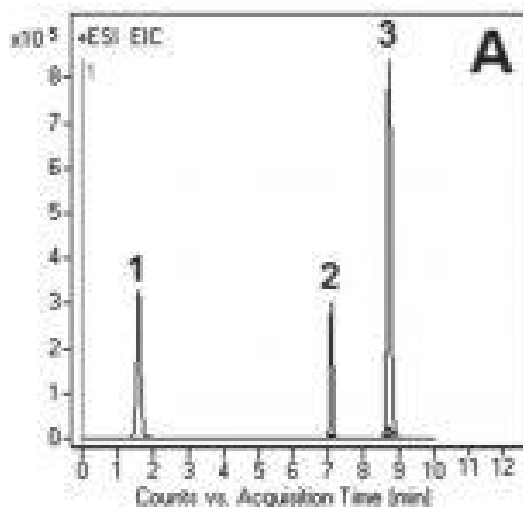

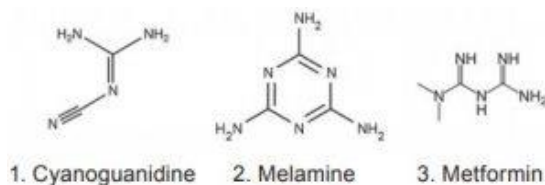

### Peaks:

1. Cyanoguanidine 85.0509 m/z (M+H)<sup>+</sup>
2. Melamine 127.0727 m/z (M+H)<sup>+</sup>
3. Metformin 130.1087 m/z (M+H)<sup>+</sup>

### Method Conditions

**Column:** Cogent Diamond Hydride™, 4μm, 100Å

**Catalog No.:** [70000-15P-2](#)

**Dimensions:** 2.1 x 150mm

**Mobile Phase:**

A: 50% Isopropanol / 50% DI Water / 0.1% Acetic Acid

B: Acetonitrile / 0.1% Acetic Acid

**Gradient:**

Time (minutes)	%B
0	100
2	100
5	20
9	20
10	100

**Post Time:** 5 minutes

**Injection vol.:** 1μL

**Flow rate:** 0.4mL / minute

**Detection:** ESI – POS - Agilent 6210 MSD TOF Mass Spectrometer

---

**Note:** *k'* of Cyanoguanidine for this Method was found to be over twice that of some Reversed Phase Methods studies that used ordinary Cyano Columns and an isocratic Mobile Phase 95% DI Water / 5% Acetonitrile. A variety of Mobile Phase additives were investigated, including 0.1% Formic Acid, 0.1% Trifluoroacetic Acid (TFA), 0.1% Trifluoroacetic Acid (TFA) +1g / L Sodium Octyl Sulfate, and 10mM Ammonium Acetate. The Reversed Phase Methods did not produce *k'* > 0.3.

---



**Attachment No 146 Impurities Determination in Metformin HCL Formulation pdf 0.3 Mb**

[Download File](#)

Printed from the Chrom Resource Center

Copyright 2025, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900

Fax: (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)