

## Method Development of Maleic Hydrazide - Tips & Suggestions

*Date: 1-OCTOBER-2012 Last Updated: 30-NOVEMBER-2025*

Maleic hydrazide has a pKa of 5.62 and a logP of -1.96. This logP value means it is very hydrophilic and will NOT retain well in Reversed Phase HPLC.

For this reason, we would suggest you try a method in ~~Aqueous Normal Phase~~ ANP using the [Cogent Diamond Hydride™](#) column in order to retain it. The pKa of 5.62 suggests that a mobile phase additive of 0.1% formic acid would work best. And 0.1% TFA can also be investigated if detection is not done by MS.



Printed from the Chrom Resource Center  
Copyright 2025, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451

Tel: (732) 380-8900  
Fax: (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)