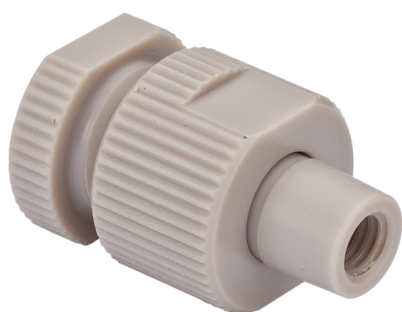


## Catalog Number

49681-TF

**Inline Filter**, for HPLC, female connections All-PEEK filter holder only. Uses disposable filters which are sold separately. ARE-Applied Research brand. Bulk Price.

**Bulk Pricing****Additional Info**

Compatible with typical flow rate analytical columns up to 4,500 psi and 5 ml/minute. Also used as a union to connect tubing, but with a filter. *Not recommended for use with UHPLC pressure ranges.*

- Economical way to protect your columns and instruments.
- Re-usable filter holder.
- Easy to install after the pump, before the injector.
- Easy to replace the filters.

## Additional Info

The universal connections (fits all standard HPLC threads and pilot lengths) can be installed with a wrench or hand tightened.

When these filters are in use, and pressure starts to rise, simply disconnect (after pressure in the column is at atmospheric), dispose of and replace with a suggested, new 0.188 inch, 0.5 um stainless steel, PEEK or Titanium PEEK-encased filter; sold separately but any filter that is 1/4 inch OD and 1/16th inch thick should work.

- Click [HERE](#) for suggested 0.5 um SS filter ordering information.
- Click [HERE](#) for economical pack of 10 PEEK holders.

### MicroSolv

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)